

# Install MediaRecorder

**IMPORTANT** - If you have an older version of MediaRecorder - Uninstall MediaRecorder 6.0 and older versions before you install MediaRecorder 6.5.

## To install MediaRecorder 6.5

1. Download the installation files from [my.noldus.com](https://my.noldus.com), and extract the contents.
2. Double-click the file **MediaRecorder [Version number] Setup.exe**.
3. Under Installation type, select **Standard**.
4. In the **Drivers and tools** field, choose the driver of your video device.
5. Click **Next**, accept the End-User License Agreement and click Install.

### Notes

- See [Specifications](#) for the supported devices.
- See [Recording Devices](#) for specific information on your device like the maximum number of cameras you can use simultaneously and the maximum supported frame rate and resolution.

---

Revision #11

Created 29 November 2023 15:23:05

Updated 20 December 2023 17:41:29