


Record video

When you selected all cameras, you see a preview in the main MediaRecorder window.

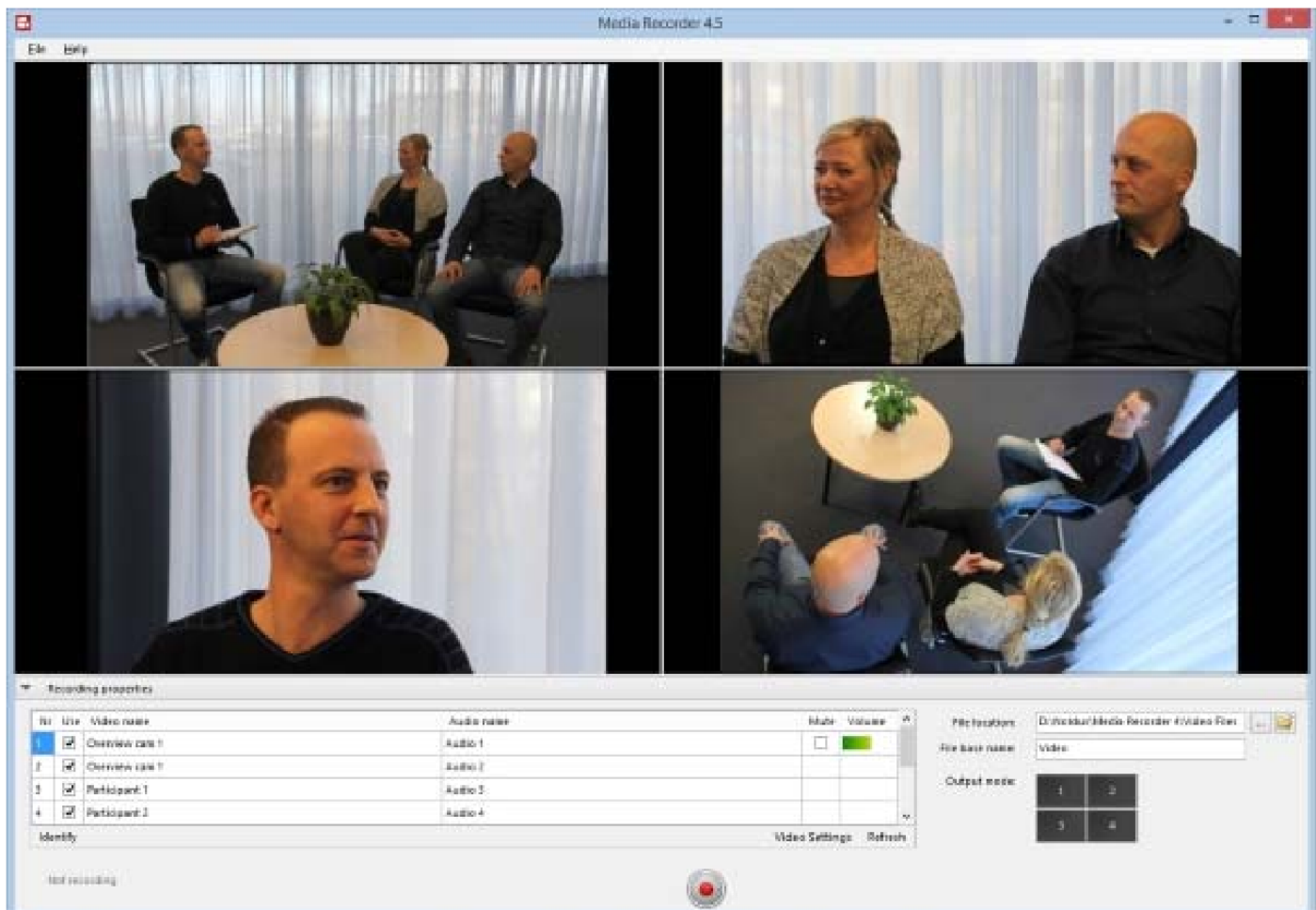

TIP If you selected multiple microphones, it can be useful to mute all but one audio channel in the preview. Select the audio source in the **Mute** column of the table under the preview. Muting audio in the preview does not influence the recording.


Nr	Use	Video name	Audio name	Mute	Volume
1	<input checked="" type="checkbox"/>	Video 1	Audio 1	<input type="checkbox"/>	<div></div>
2	<input checked="" type="checkbox"/>	Video 2	Audio 2	<input checked="" type="checkbox"/>	<div></div>
3	<input checked="" type="checkbox"/>	Video 3	Audio 3	<input checked="" type="checkbox"/>	<div></div>
4	<input type="checkbox"/>	Video 4	Audio 4	<input type="checkbox"/>	

IdentifyVideo SettingsRefresh

To start recording, press the **Start recording** button.

	Mute	Volume
	<input checked="" type="checkbox"/>	

Video Settings Refresh

A circular button with a red center and a grey outer ring, highlighted by an orange square border.

A pop-up window will appear, informing you about Noldus's Data Transparency Policy. Read it carefully before clicking the **I will inform the participants** button. You can also check the box **Don't show me again** to prevent further occurrence.

Data transparency

Responsible Use

Informed consent

You are starting the recording. Before you continue, if your study involves human subjects, please inform the participants in your study what you will be doing with their data and why. Make sure you have their consent to take part in the research.

Data Transparency

By clicking the OK button, you acknowledge your responsibility to inform the participants about:

- Which data is collected
- How the data will be used
- Where the data will be stored
- Who will have access to the data
- What the collected data will not be used for
- How the participant will benefit from the data collection

Use of MediaRecorder is subject to certain restrictions, which are documented in section 7 of the [End-User License Agreement](#).

☐ Don't show me again


I will inform the participants

The button then changes into the **Stop Recording** button and the Recording timer starts running.

2	<input checked="" type="checkbox"/>	Participant 1	Audio 3
3	<input checked="" type="checkbox"/>	Overview cam 1	Audio 2
4	<input checked="" type="checkbox"/>	Participant 2	Audio 4

Identify


Recording: 00:00:03 (hh:mm:ss)



Press the button again to stop recording. Recording stops automatically when you [specified a recording time](#) .

how to access the recorded videos


To quickly access the recorded videos, press the **Open target folder** button in the main MediaRecorder window.

File location: ... 

File base name:

Output mode:

1	2
3	



The video files are stored in the default video file location, or the location you specified. See [To select the video file location](#).

Revision #8

Created 29 November 2023 15:49:07

Updated 11 March 2024 10:01:18