

Carry out an observation in The Observer XT

Procedure

1. In The Observer XT, choose **Observe > Observation > New** to create an observation.
2. See **Carry out your observation** in The Observer XT Help.
3. Press the **Start observation** button and score the data. See **Start the observation** and **Score data** in The Observer XT Help.
4. Press the **Stop observation** button. See **Stop the observation** in The Observer XT Help.

Note

IMPORTANT Allow enough time between the actions. If you start an observation before MediaRecorder is completely open, video recording will not start.

Revision #4

Created 29 November 2023 22:23:38

Updated 20 December 2023 14:29:04 by Viktor Kovacs