

# Carry out an observation in The Observer XT

## Procedure

1. In The Observer XT, choose **Observe > Observation > New** to create an observation.
2. See **Carry out your observation** in The Observer XT Help.
3. Press the **Start observation** button and score the data. See **Start the observation** and **Score data** in The Observer XT Help.
4. Press the **Stop observation** button. See **Stop the observation** in The Observer XT Help.

## Note

**IMPORTANT** Allow enough time between the actions. If you start an observation before MediaRecorder is completely open, video recording will not start.

---

Revision #4

Created 29 November 2023 23:23:38

Updated 20 December 2023 15:29:04 by Viktor Kovacs