

# Change the video file location

## Aim

To save the videos in another location than the default one.

The default location where MediaRecorder stores its video files is:

D:\Noldus\MediaRecorder\Video Files — If Noldus IT installed your system


C:\Users\Public\Public Documents\Noldus\MediaRecorder\Video Files — In all other cases. It is recommended to set this folder to be located on a separate, fast, large capacity drive.

## Prerequisite

You must be able to write in the selected folder and there must be enough file space available (at least 600 Mb per video per hour).

## Procedure

Click this button and browse to the new location.

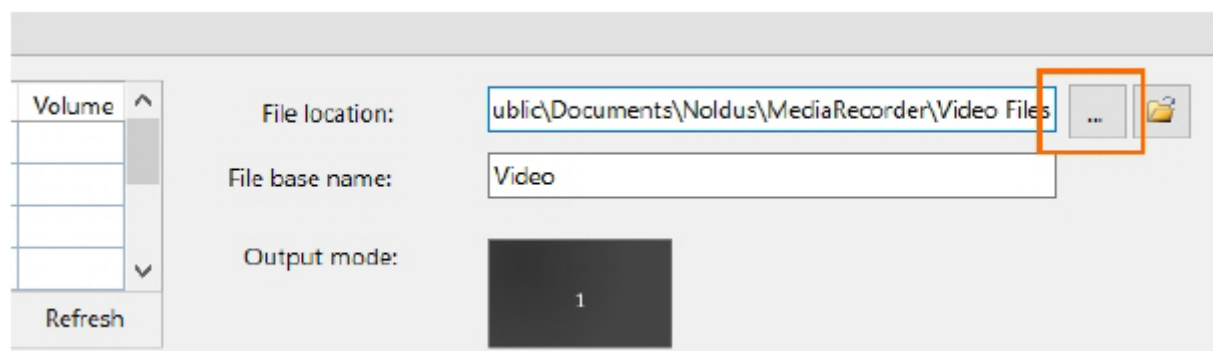

**tip** Click this button to open the folder with video files in Windows Explorer and play recorded videos.

