

# Create audio settings

## Aim


To optimize the recording volume.

## Prerequisites

- You have Axis IP cameras that are supported with MediaRecorder.  
See [Supported IP cameras](#) in [Recording devices](#)
- You followed the procedure in [Ethernet card setup](#)
- You connected all cameras.
- You gave the cameras an IP address.  
See [Set camera IP address](#)
- You gave the camera a password and selected the correct power line frequency.  
See [Set camera password and select power line frequency](#)
- You followed the procedure in [Create camera settings](#)

## Procedure

1. Open the camera settings,  
See [Access camera settings](#)
2. Open the **Audio** tab.
3. Adjust the **Gain** slider and check in the **Input** section that the green field in the **Level** field is maximal with normal speech.

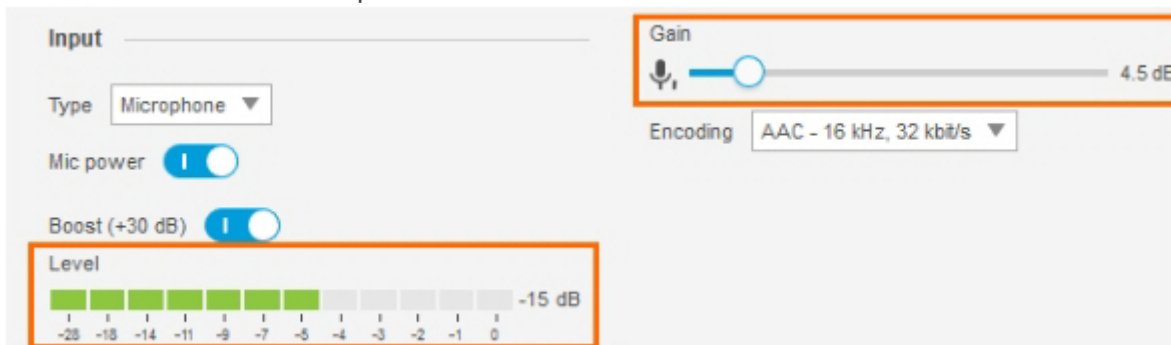

## Note

You cannot record the audio of non-ONVIF devices in MediaRecorder. Optionally record audio from a separate microphone that is connected to the microphone input of the computer with MediaRecorder. The Axis cameras supported with MediaRecorder support ONVIF.

Revision #6

Created 1 December 2023 09:49:17

Updated 20 December 2023 13:47:08