

# Create MediaRecorder Settings

## Aim

To connect MediaRecorder to N-Linx Server and select the video file folder.

## Prerequisites

- If MediaRecorder, The Observer XT, and N-Linx Server are not all installed on the same computer, you created exceptions for the N-Linx port Windows Firewall.


See [Create exceptions for N-Linx port in Windows Firewall](#)

- You installed N-Linx Server on a computer in the network.
- N-Linx Agent is installed on the computer with MediaRecorder and connected to N-Linx Server.

See [Install N-Linx Agent](#)

## Procedure

1. In MediaRecorder, choose **File > Preferences > N-Linx settings**

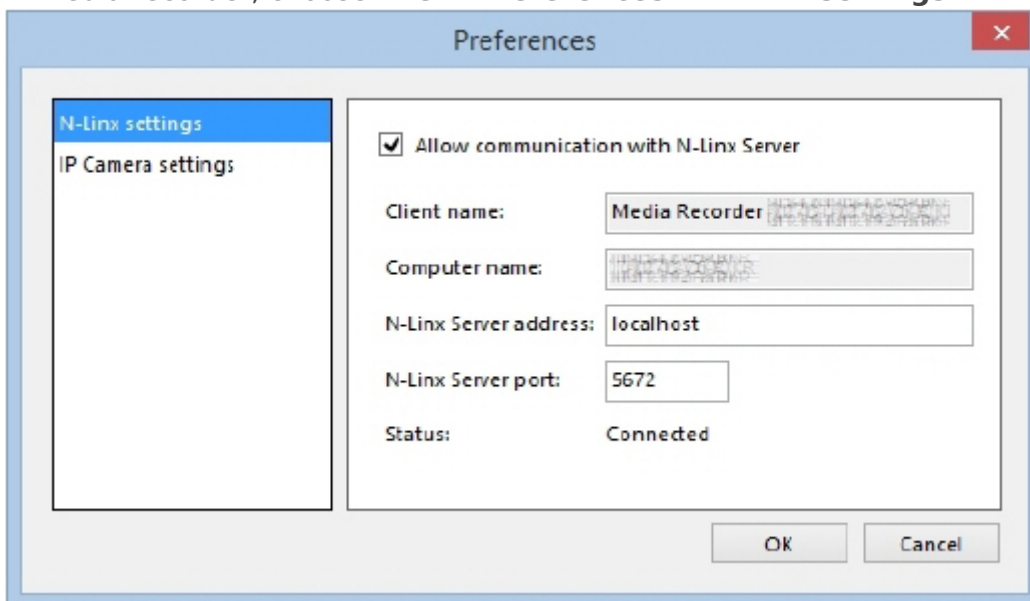

2. Select **Allow communication with N-Linx Server**.
3. In the **N-Linx Server address** field, leave the default value Localhost if MediaRecorder runs on the same computer as N-Linx Server.  
If the programs run on different computers, enter the name of the computer with N-Linx Server. See also [To find the computer name](#)
4. In the **N-Linx Server port** field, leave the default port 5672. Ask your system administrator for assistance if this port is used by another program.

5. When the connection with N-Linx Server is established the **Status** will be **Connected**. Click **OK**.
6. If MediaRecorder and The Observer XT run on different computers, select the mapped drive as location to store the video files and not the original folder. The original folder is locally on the computer with MediaRecorder and cannot be accessed by The Observer XT.

**example** You mapped the default folder C:\Users\Public Documents\Noldus\MediaRecorder\Video files as **X:\**. Set the default video file location where MediaRecorder stores its video files to **X:\** and not to the original folder.

---

Revision #6

Created 29 November 2023 23:19:15

Updated 20 December 2023 13:47:09