


Enter a default user name and password of IP cameras

Aim

To enter the ONVIF user name and ONVIF password of IP cameras only once, in case you have several cameras with the same credentials.

Procedure

1. Choose **File > Preferences > IP camera settings**.
2. Select **Use default IP camera user name and password**.
3. Enter the user name and password.

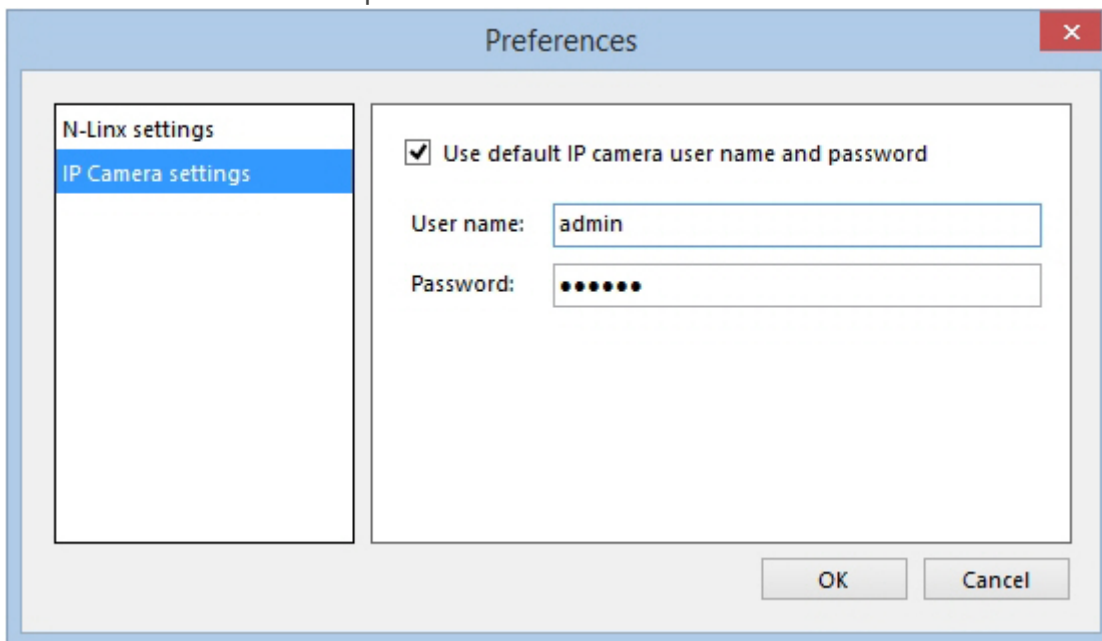

4. Follow the procedure in [Select ONVIF cameras in MEdiaRecorder](#) or [Select devices that do not support ONVIF in MediaRecorder](#). When you click the **Advanced video settings** button, the user name and password are filled in automatically.

Revision #5

Created 29 November 2023 22:50:45

Updated 20 December 2023 13:47:09