


# Enter a default user name and password of IP cameras

## Aim

To enter the ONVIF user name and ONVIF password of IP cameras only once, in case you have several cameras with the same credentials.

## Procedure

1. Choose **File > Preferences > IP camera settings**.
2. Select **Use default IP camera user name and password**.
3. Enter the user name and password.

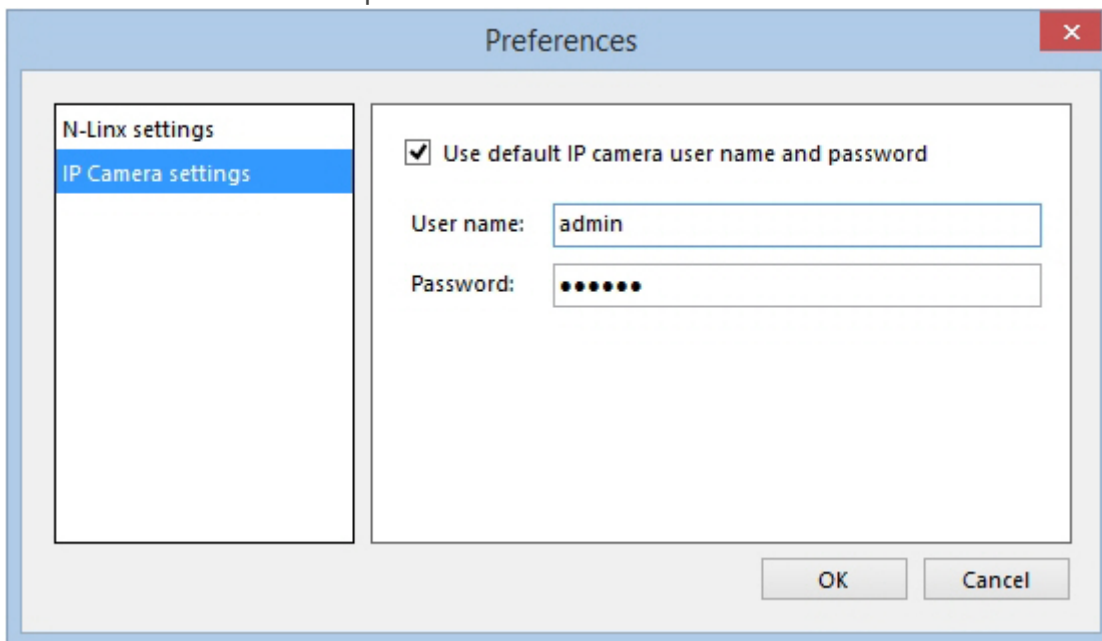

4. Follow the procedure in [Select ONVIF cameras in MEdiaRecorder](#) or [Select devices that do not support ONVIF in MediaRecorder](#). When you click the **Advanced video settings** button, the user name and password are filled in automatically.

---

Revision #5

Created 29 November 2023 21:50:45

Updated 20 December 2023 12:47:09