

# Install MediaRecorder and device drivers

## Aim


To install MediaRecorder and, for some recording devices, their drivers.

## Procedure

**IMPORTANT** - If you have an older version of **MediaRecorder** - Uninstall MediaRecorder 6.0 and older versions before you install MediaRecorder 6.5.

1. Download the installation files from [my.noldus.com](http://my.noldus.com), and extract the contents.
2. Browse to the file **MediaRecorder [Version number] Setup.exe** and double-click it.

3. We recommend to choose **Standard** as Installation type. Choose **Custom** only to change the location where the program is installed.

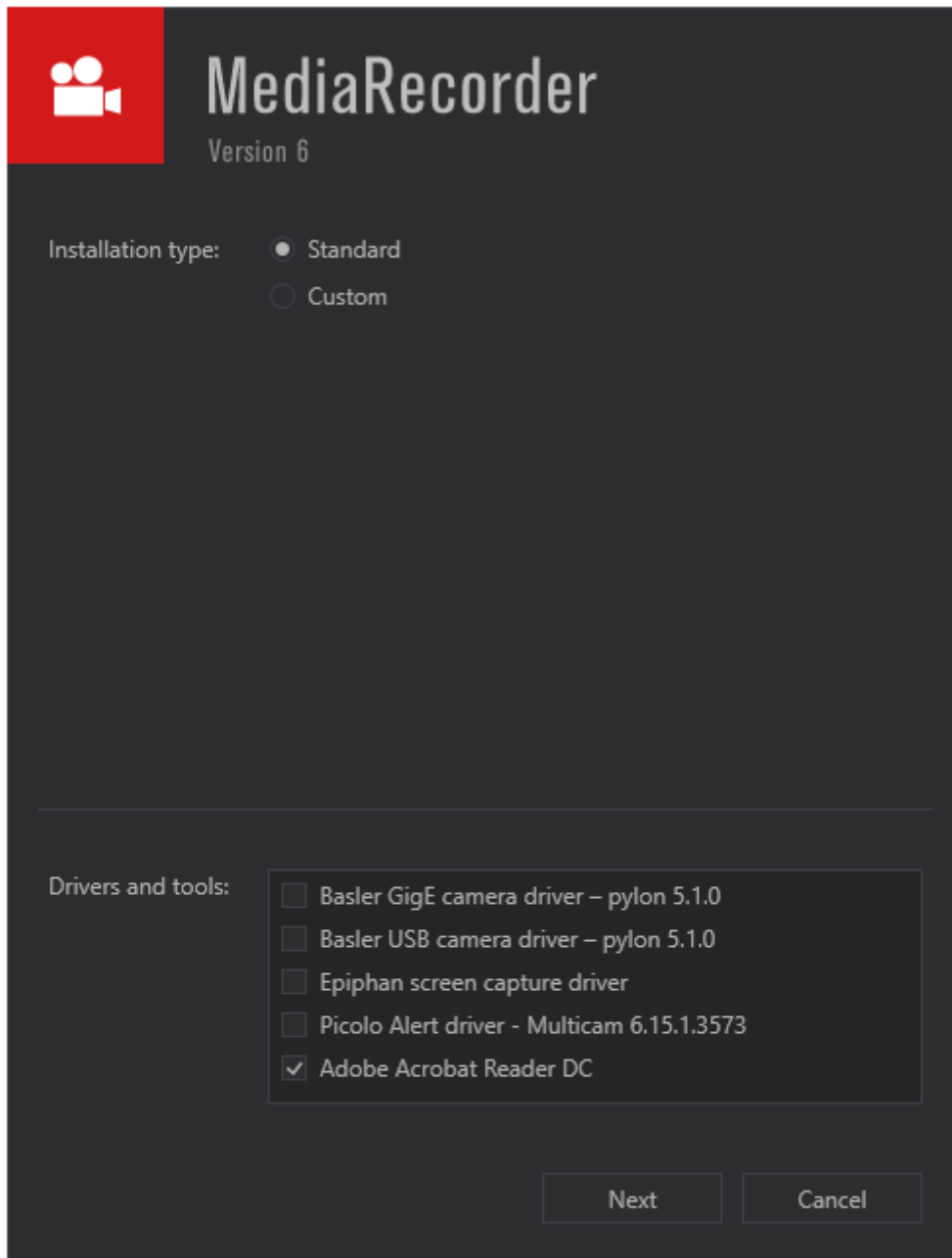

4. In the **Drivers and tools** field, choose the driver of your video device.
5. Click **Next**, accept the License Agreement and click **Install**.

The following setup prerequisites are installed automatically before MediaRecorder is installed:

- Sentinel Run time — Needed for correct functioning of your hardware key.
- Noldus LeadTools Package 21.
- Noldus MediaLooks Proxy Filters 2.1 — Needed for screen capture of the monitor of the computer with MediaRecorder.
- Medialooks Screen Capture 2.0.3 — Needed for screen capture of the monitor of the computer with MediaRecorder.
- Microsoft .NET Framework 4.7.1 —To run MediaRecorder. This program is installed only when not already present on the computer.

Revision #6

Created 29 November 2023 17:12:43

Updated 20 December 2023 22:26:42