

Install the drivers for the Picolo Alert card

Aim

To install the software that allows communication between the Euresys card and the computer operating systems.

Prerequisites

- You use MediaRecorder with analog cameras.
- You did not buy a complete setup from Noldus IT.
- You installed the Euresys Picolo Alert card into the computer.

Procedure

1. Insert the MediaRecorder installation USB stick into your computer.
2. If you have the Picolo Alert card, open the folder Drivers and Tools\Drivers\Euresys\Alert and double-click the file **multicam-win10-6.15.1.3573.exe**
3. Follow the instructions on your screen.
4. When installation is finished, a message appears that the drivers were successfully installed. Click **Close**.
5. Installing the Picolo card drivers involves a a computer restart. After restart, the installation is finished.

Continue with installing MediaRecorder. See [Installation](#)

Revision #5

Created 29 November 2023 23:50:53

Updated 20 December 2023 13:47:08