

Mute audio in preview

Aim

To mute audio sources in MediaRecorder. The muted audio is recorded normally in the video files.

Prerequisite

You selected microphones in the **Video Settings** window.

See [Select an audio device](#)

Procedure

Select the audio sources you wish to mute in the **Mute** column.

| Nr | Use | Video name | Audio name | Mute | Volume |
|---------------------------------|-------------------------------------|------------|------------|-------------------------------------|-------------|
| 1 | <input checked="" type="checkbox"/> | Video 1 | Audio 1 | <input type="checkbox"/> | <div></div> |
| 2 | <input checked="" type="checkbox"/> | Video 2 | Audio 2 | <input checked="" type="checkbox"/> | <div></div> |
| 3 | <input checked="" type="checkbox"/> | Video 3 | Audio 3 | <input checked="" type="checkbox"/> | <div></div> |
| 4 | <input type="checkbox"/> | Video 4 | Audio 4 | <input type="checkbox"/> | <div></div> |
| Identify Video Settings Refresh | | | | | |

Revision #6

Created 29 November 2023 22:39:26

Updated 20 December 2023 23:59:55