

# Mute audio in preview

## Aim

To mute audio sources in MediaRecorder. The muted audio is recorded normally in the video files.


## Prerequisite

You selected microphones in the **Video Settings** window.

See [Select an audio device](#)

## Procedure

Select the audio sources you wish to mute in the **Mute** column.

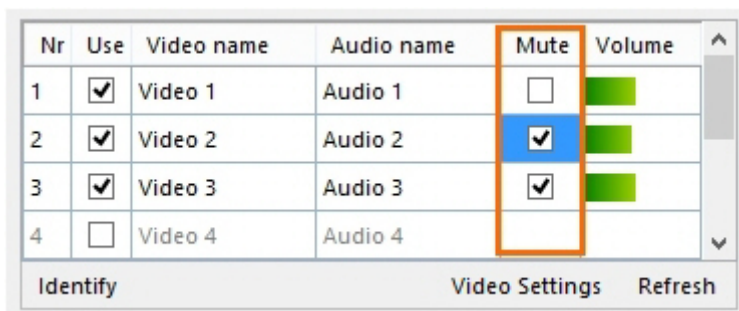

Nr	Use	Video name	Audio name	Mute	Volume
1	<input checked="" type="checkbox"/>	Video 1	Audio 1	<input type="checkbox"/>	<div></div>
2	<input checked="" type="checkbox"/>	Video 2	Audio 2	<input checked="" type="checkbox"/>	<div></div>
3	<input checked="" type="checkbox"/>	Video 3	Audio 3	<input checked="" type="checkbox"/>	<div></div>
4	<input type="checkbox"/>	Video 4	Audio 4	<input type="checkbox"/>	<div></div>

Identify Video Settings Refresh

Revision #6

Created 29 November 2023 22:39:26

Updated 20 December 2023 23:59:55