

Record your videos

Aim


To create video files from the video stream of the connected recording devices.

Prerequisites


- Your devices are connected to the computer with MediaRecorder, and, if applicable, switched on.
- MediaRecorder is open.
- You selected the cameras and, optionally, microphones and specified the output of the video files in the **Video Settings** window (**File > Video Settings**).

Procedure


1. When you selected all settings in the **Video Settings** window, click **OK** to return to the Main window. You will see a preview of the cameras.

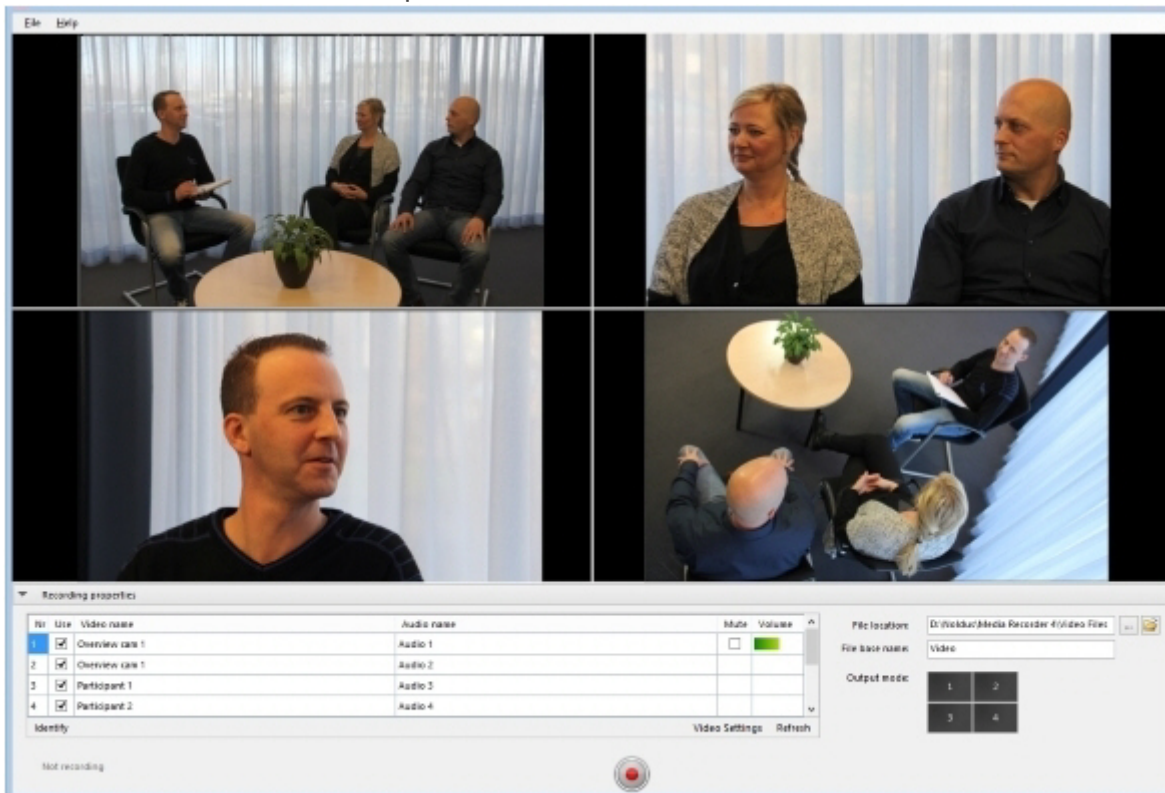


2. Optionally, adjust options like the video file location, the video file base name or the order of the video files. See [Options](#)
3. Press the **Start recording** button.

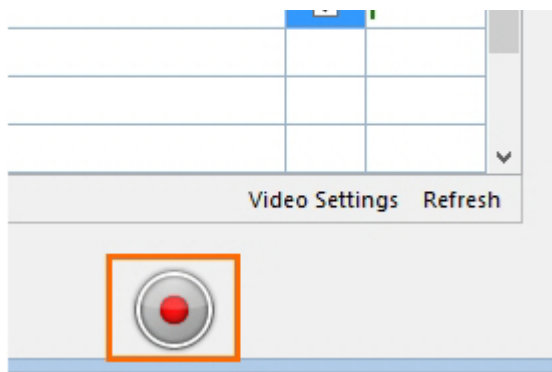

- A pop-up window will appear, informing you about Noldus's Data Transparency Policy. Read it carefully before clicking the **I will inform the participants** button. You can also check the box **Don't show me again** to prevent further occurrence.

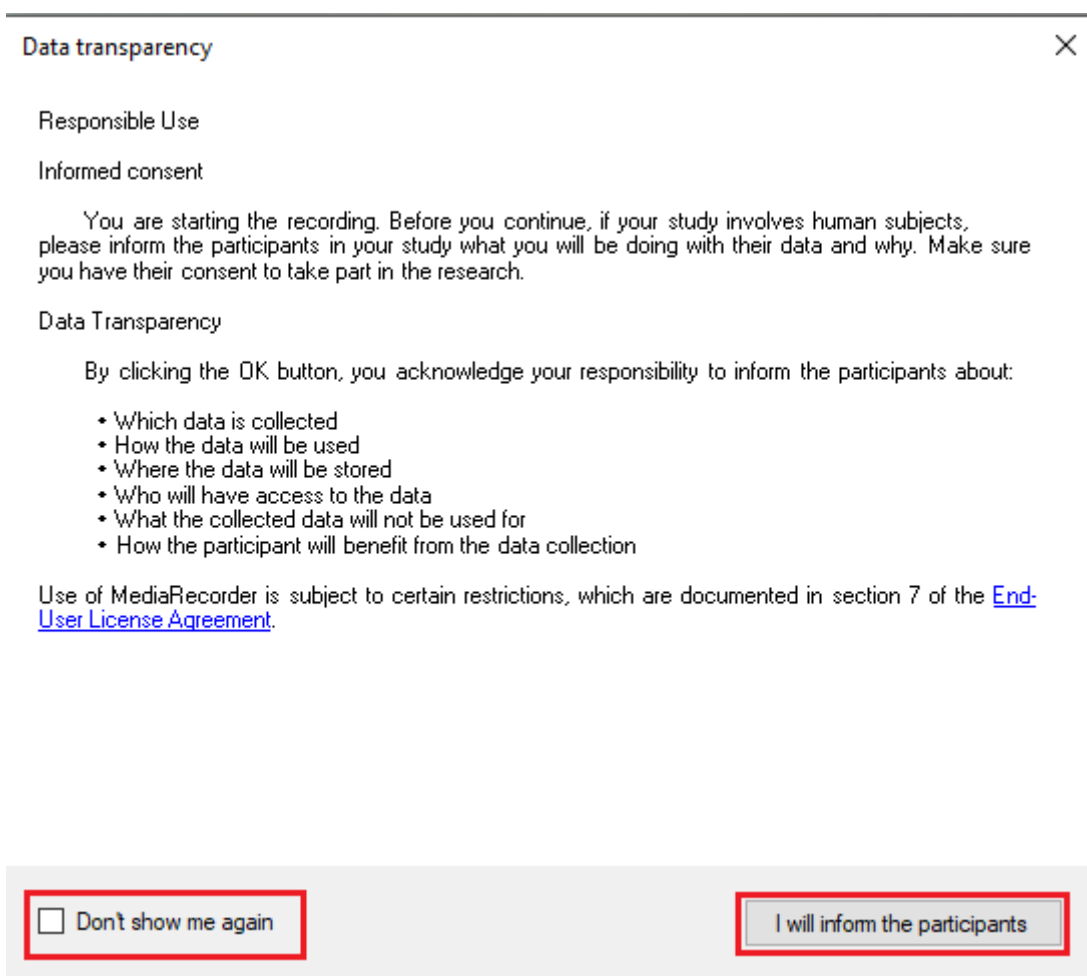

- The button changes into the **Stop Recording** button and the Recording timer starts running.

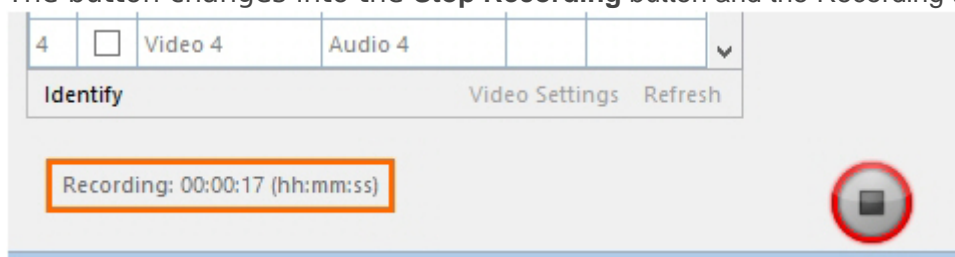

6. Press the button again to stop recording. Recording stops automatically when you specified a [Maximum recording time](#).

Note

If you have IP cameras, see [Pan, Tilt, Zoom control with IP cameras](#) how to change the camera view before or during recording.

Revision #9

Created 29 November 2023 22:34:26

Updated 20 December 2023 23:58:47