

# Record your videos

## Aim


To create video files from the video stream of the connected recording devices.

## Prerequisites


- Your devices are connected to the computer with MediaRecorder, and, if applicable, switched on.
- MediaRecorder is open.
- You selected the cameras and, optionally, microphones and specified the output of the video files in the **Video Settings** window (**File > Video Settings**).

## Procedure

1. When you selected all settings in the **Video Settings** window, click **OK** to return to the Main window. You will see a preview of the cameras.

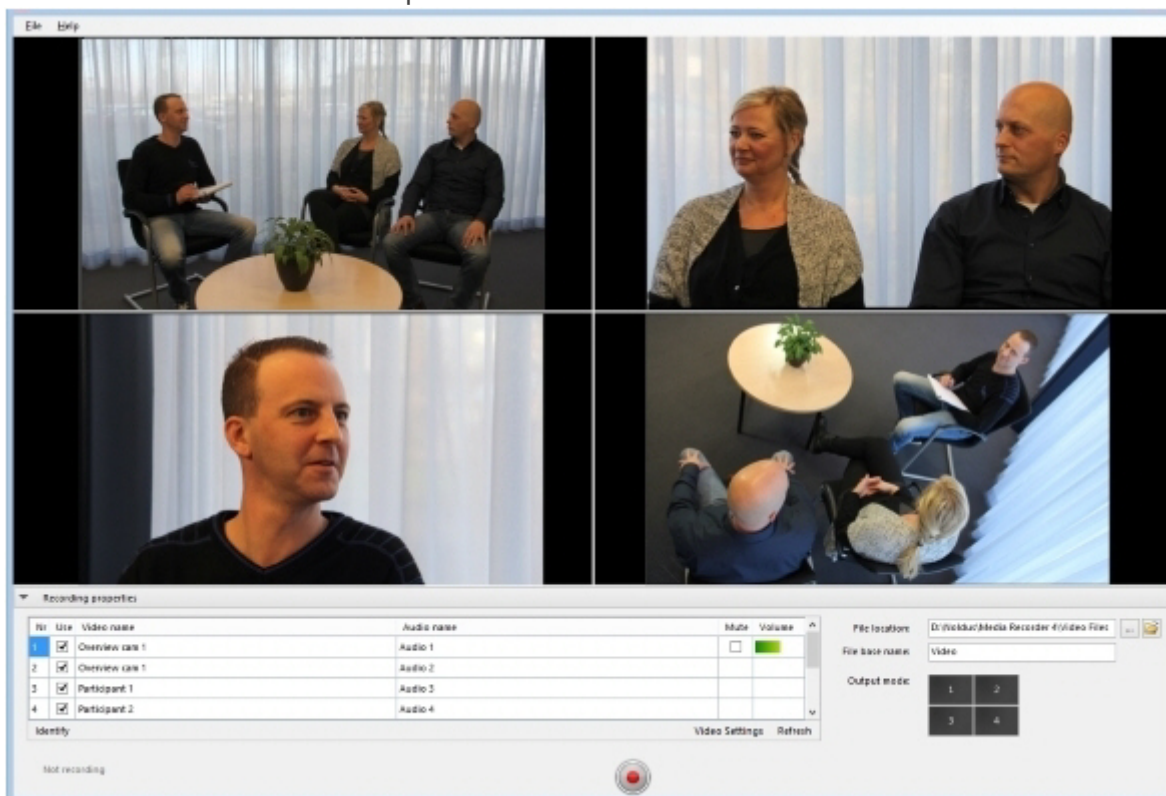

2. Optionally, adjust options like the video file location, the video file base name or the order of the video files. See [Options](#)
3. Press the **Start recording** button.

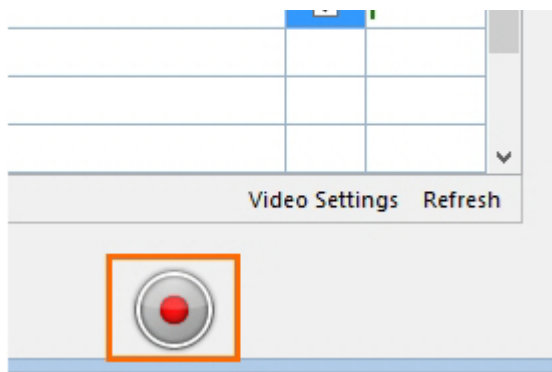

- A pop-up window will appear, informing you about Noldus's Data Transparency Policy. Read it carefully before clicking the **I will inform the participants** button. You can also check the box **Don't show me again** to prevent further occurrence.

Data transparency

Responsible Use

Informed consent

You are starting the recording. Before you continue, if your study involves human subjects, please inform the participants in your study what you will be doing with their data and why. Make sure you have their consent to take part in the research.

Data Transparency

By clicking the OK button, you acknowledge your responsibility to inform the participants about:


- Which data is collected
- How the data will be used
- Where the data will be stored
- Who will have access to the data
- What the collected data will not be used for
- How the participant will benefit from the data collection

Use of MediaRecorder is subject to certain restrictions, which are documented in section 7 of the [End-User License Agreement](#).

☐ Don't show me again

I will inform the participants

- The button changes into the **Stop Recording** button and the Recording timer starts running.

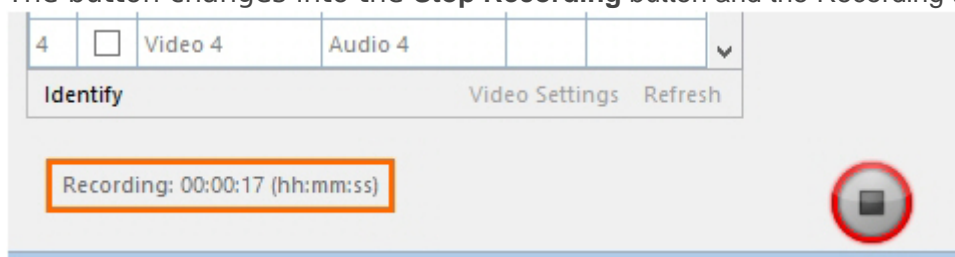

6. Press the button again to stop recording. Recording stops automatically when you specified a [Maximum recording time](#).

## Note

If you have IP cameras, see [Pan, Tilt, Zoom control with IP cameras](#) how to change the camera view before or during recording.

---

Revision #9

Created 29 November 2023 21:34:26

Updated 20 December 2023 22:58:47