

# Set video file base name

## Aim

To select a word or phrase to start the video file name with.

The video file name consists of the **File base** name followed by:


- The video input name.
- The date and the time the videos were recorded.
- The number of the video camera.

Example – The **File base name** is Video, you changed the name of Video 2 to Webcam in the **Video Settings** window (**File > Video Settings**), the date is 11 May 2016, the time 10:54:43 and you create 2 videos simultaneously. The name of the second video will then be:

Video Webcam 5\_11\_2016 10\_54\_43 AM 2.mp4

## Procedure

Enter a name in the **File base name** field.

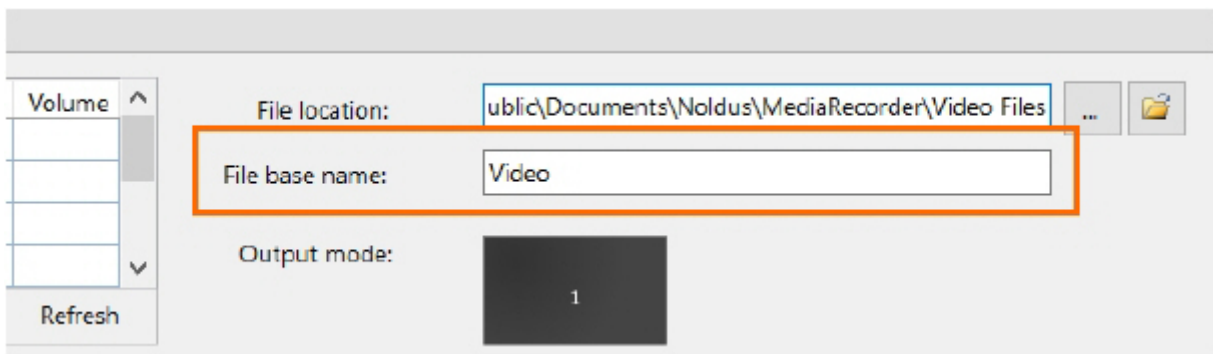

The screenshot shows a software window titled 'Video Settings'. On the left is a 'Volume' control with a slider and a 'Refresh' button. The main area contains three settings: 'File location' with a text box showing 'ublic\Documents\Noldus\MediaRecorder\Video Files' and a folder icon button; 'File base name' with a text box containing 'Video', which is highlighted by an orange rectangular box; and 'Output mode' with a dark grey button labeled '1'.

## Note

The date and time depend on your computer's regional settings. To change these, open the **Control panel** and then **Region and Language**. Select your preferences in one of the tabs.

---

Revision #3

Created 29 November 2023 22:37:59

Updated 21 December 2023 00:01:20