


Show frame rate, resolution and format

Aim

To display hidden rows in the table with video devices.

Procedure

Right-click one of the column headers and select the preferred options.

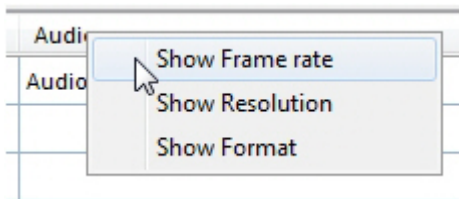

Notes

- The **Format** column shows the color space of the video file. The column is hidden by default, because the optimal color space is selected by default. See the following page for more information on color space.
<http://www.noldus.com/knowledge-base/what-color-space-format-used-media-recorder>.
- **TIP** Follow the same procedure to show or hide the Frame rate, Resolution, and Format columns in the **Video Settings** window.

Revision #3

Created 29 November 2023 22:40:33

Updated 21 December 2023 00:10:40