

Upgrade MediaRecorder

Aim

To upgrade an older version of the MediaRecorder software to version 6.5.

Prerequisites

You have an older MediaRecorder version.

You obtained an activation key for your MediaRecorder license.

Procedure

1. Download the installation files from my.noldus.com, and extract the contents.
2. Browse to the file **MediaRecorder [Version number] Setup.exe** and double-click it.
3. Follow the steps on your screen to install MediaRecorder 6.5.
4. [Activate your MediaRecorder license](#).

Revision #9

Created 29 November 2023 17:07:09

Updated 21 December 2023 10:20:34 by Niek Wilmink