

# Use of the Screen Capture tool

## Using the standard Screen Capture tool in MediaRecorder

The Screen Capture tool is an add-in to your MediaRecorder license. It allows you to record the screen of the computer with MediaRecorder. It is a software tool that is built in in MediaRecorder.

**IMPORTANT** To record the screen of a computer other than the one with MediaRecorder installed, use an Epiphan screen capture device instead.

Prerequisites

- You record the full screen of the PC with MediaRecorder.
- You record one screen only.
- You use MediaRecorder on the Dell Precision 3551 mobile workstation.


See also [System requirements](#).

## Drivers

For the MR screen capture option, you do not need to install device drivers.

## Supported setup (portable lab)

The table below shows the maximum supported (1) resolution, (2) frame rate (fps), and (3) recording time (hours) on a 3551 notebook.

	Device	Resolution	Frame rate	Recording time
	Logitech Brio + Built-in screen capture	1920 x 1080	30	1

Other USB cameras may also work. However, it could be that you have to change the Color space to MJPG. See [Show frame rate, resolution and format](#).

You can also combine built-in screen capture with, for instance, an IP camera.

## Procedure


1. In MediaRecorder, choose **File > Video Settings**.
2. From the **Video device** list, select **Medialooks Screen Capture**.

	Use	Video name	Video device		Frame rate	Resolution
1	<input checked="" type="checkbox"/>	Video 1	Medialooks Screen Captu		30.00	720 x 486
2	<input type="checkbox"/>	Video 2	USB Video Device			
3	<input type="checkbox"/>	Video 3	RTSP IP Camera (Noldus RTSP Source F			
4	<input type="checkbox"/>	Video 4	Medialooks Screen Capture			
			Select video device...			

3. Click the **Advanced Video Settings** button to select recording options.

Settings						
	Use	Video name	Video device		Frame rate	Resolution
1	<input checked="" type="checkbox"/>	Video 1	Medialooks Screen Capture		30.00	1280 x 800
2	<input type="checkbox"/>	Video 2	Select video device...			

4. In the **Medialooks Screen Capture Properties** window that opens, select the frame rate and resolution in the **Output** field. Although you are allowed to select a high frame rate, the actual frame rate of the recording will depend on the processor speed of your computer. Most likely this will be in the order of magnitude of 10 frames per second.

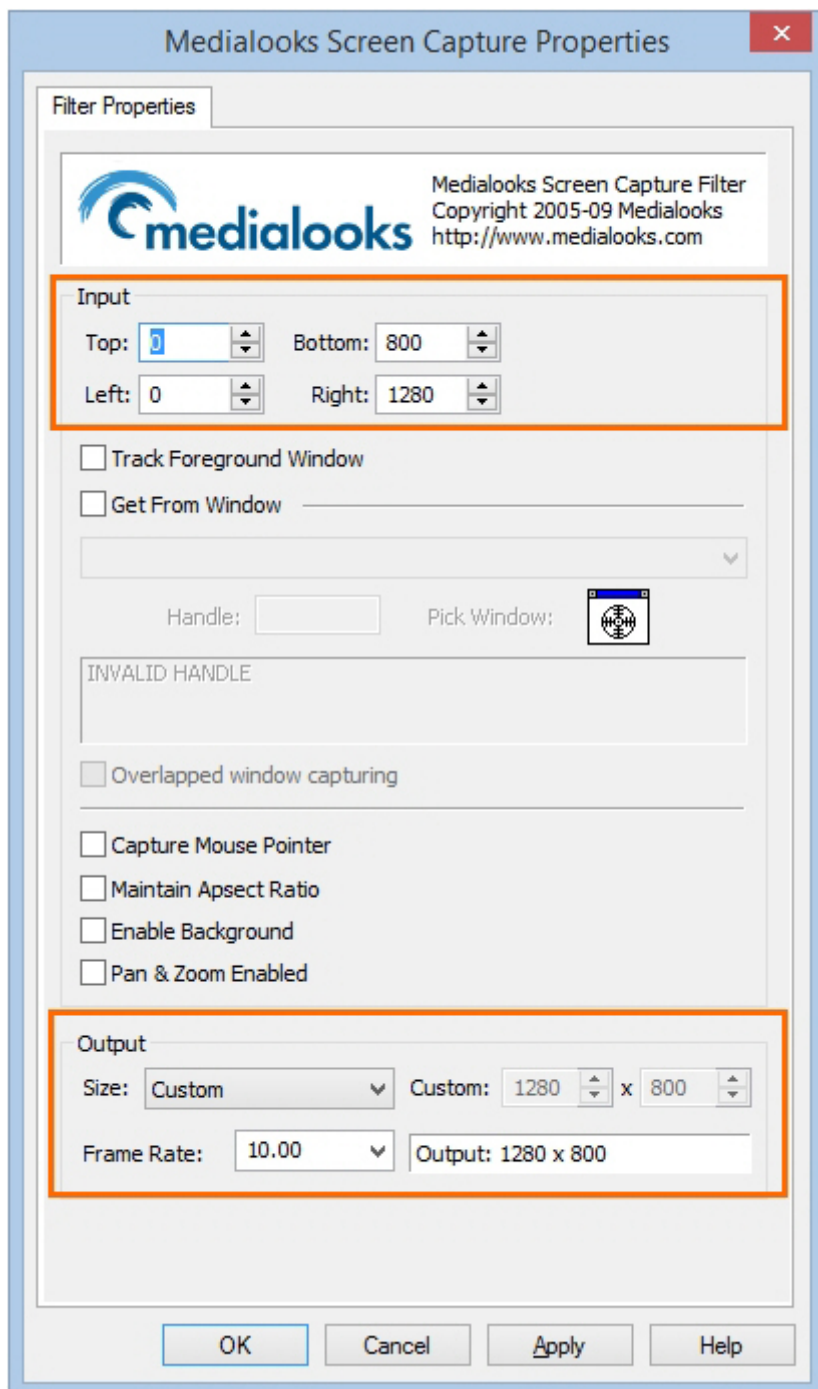

5. Although this window has several recording options, the use of the built-in screen capture is designed and supported to record one window on the MediaRecorder computer only. Minimize MediaRecorder after you start recording.

## To re-enable settings

As soon as you select the screen capture device, it is used for the video preview and you cannot change its settings. The Output fields are grayed out. To enable these fields again:

1. Close the **Medialooks Screen Capture Properties** window.
2. From the **Video device list** choose **Select video device** and click **Apply**.

	Use	Video name	Video device	Frame rate
1	<input checked="" type="checkbox"/>	Video 1	Medialooks Screen Captu ▼	30.00
2	<input type="checkbox"/>	Video 2	USB Video Device	30.00
3	<input type="checkbox"/>	Video 3	RTSP IP Camera (Noldus RTSP Source Fi	
			Medialooks Screen Capture	
4	<input type="checkbox"/>	Video 4	Select video device...	
			Select video device...	

Reselect the **Medialooks Screen Capture** device again and click the video icon to re-open the Medialooks Screen Capture Properties window (see step 3 above). You can now change the frame rate and resolution again.

## Note

The actual resolution of the recorded file may be lower than the value set in Output.

The use of built-in screen capture option is supported with a Windows text size of 100% or 125% (Control Panel > Display).

### Change the size of all items

Make text and other items on the desktop smaller and larger. To temporarily enlarge just part of the screen, use the [Magnifier](#) tool.


- ☒ **S**maller - 100%
- ☐ **M**edium - 125%
- ☐ **L**arger - 150%



### Custom sizing options

If you change Windows text size to 125%, this may result in MediaRecorder capturing only part of the screen. To solve this, do the following:

1. Make sure MediaRecorder is not running.
2. Open the folder **C:\Program Files\Noldus\MediaRecorder 6**.
3. Right-click the file **MediaRecorder.exe** and select **Properties**.
4. Open the **Compatibility** tab.
5. Select **Disable display scaling on high DPI settings** and click **Apply**.

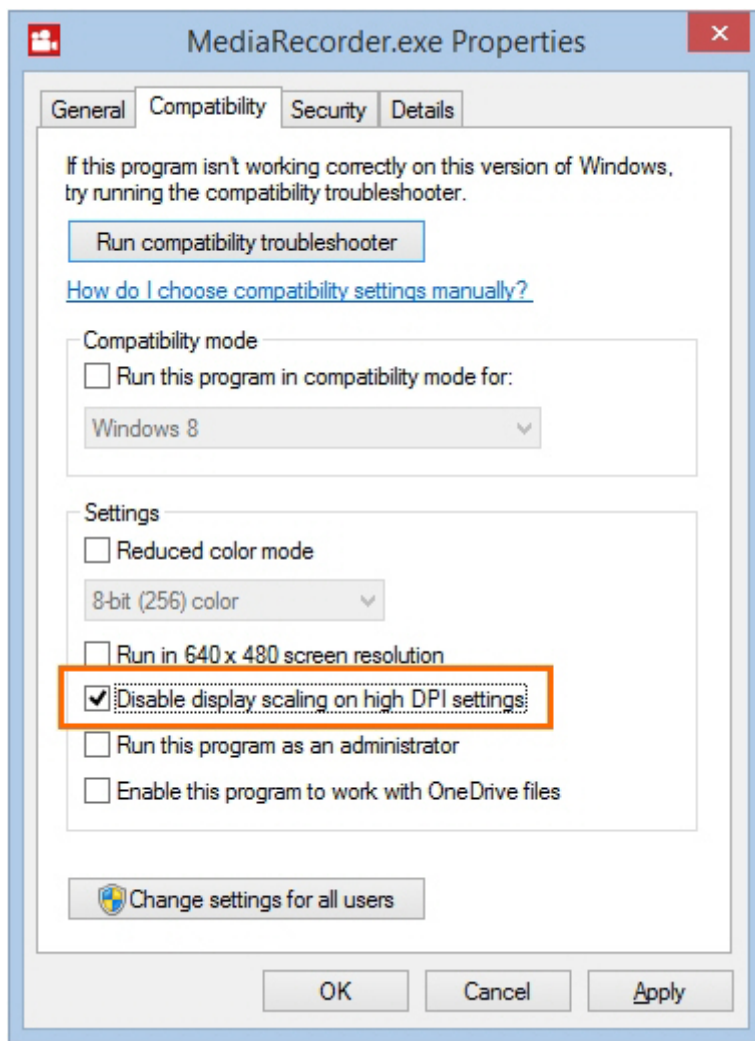

Revision #11

Created 29 November 2023 18:36:07

Updated 21 December 2023 10:25:27 by Niek Wilmink