

# View the automatic offset of videos imported into The Observer XT


## Procedure

The procedure depends on what you did:

- [You controlled video recording with The Observer XT](#)
- [You manually started video recording within 10 seconds after starting the observation in The Observer XT.](#)


## You controlled video recording with The Observer XT

Right-click the video in The Observer XT and choose **Properties**. The field **Recording details** shows the offset.

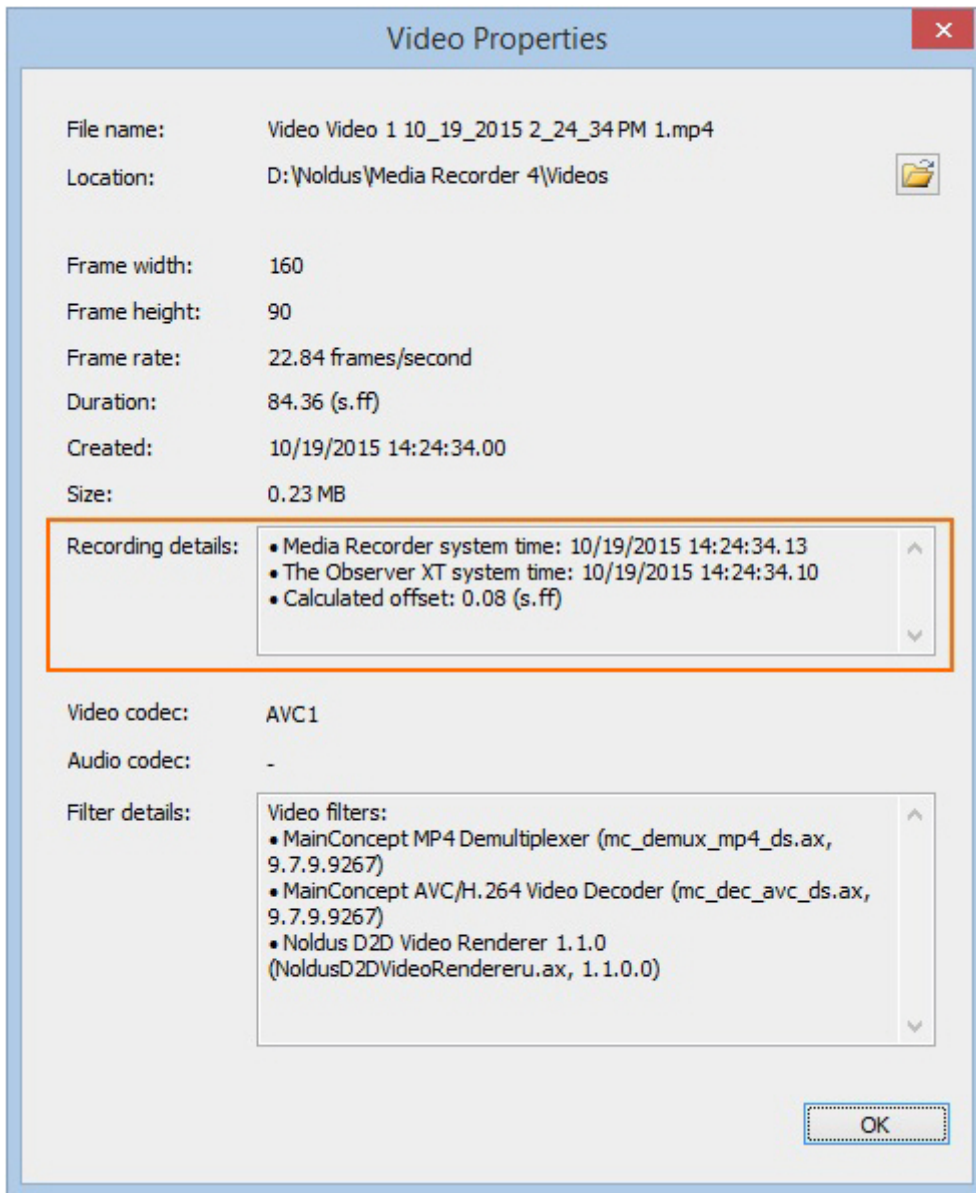


## You manually started video recording within 10 seconds after starting the observation in The Observer XT

When you import the videos into The Observer XT, an offset is automatically set. A message with the calculated offset is shown.

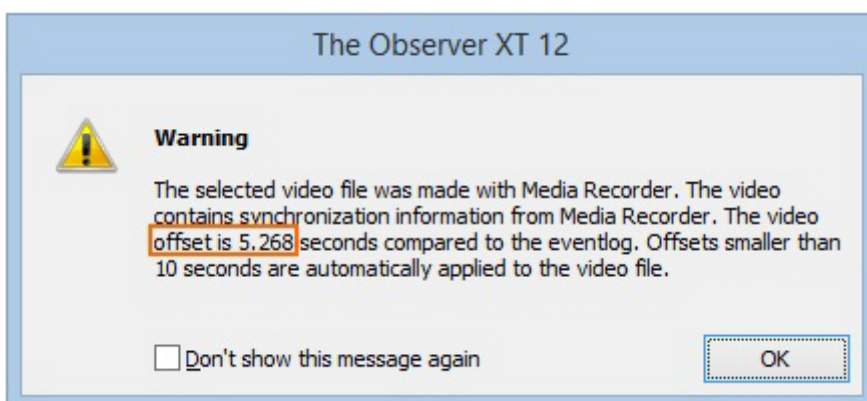

If you right-click the video in The Observer XT, the offset is also shown next to **Recording details** in the **Video Properties** window that opens.

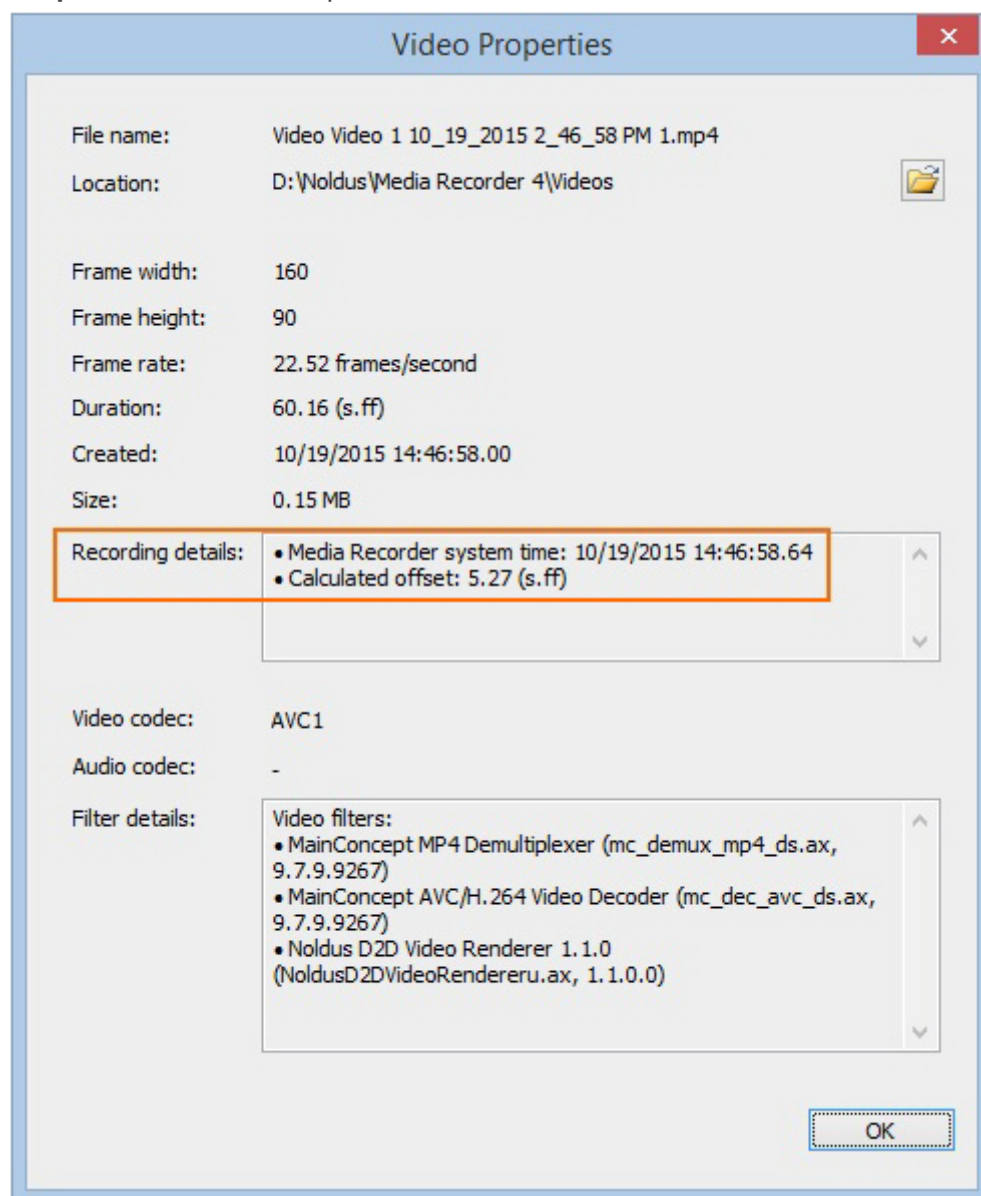

However, there can still be up to 1 second asynchronism between the observation and the video files. If that is a problem for the accuracy of your experiments, [Synchronize manually](#) to fine-tune the synchronization.

Revision #5

Created 29 November 2023 23:25:44

Updated 20 December 2023 13:47:09