

# Enter a remark

## aim

To mark something of interest that you did not define in the Codes list. This can be done during session recording or afterwards while playing the session back.

## prerequisites

- Your license includes the Annotation Module.
- To add remarks during session recording, your user role allows recording (Technician, Administrator, or Trainer).
- To add a remark while playing a session back, you recorded the session, received ownership or annotator rights for the session.

## Remarks in a multiple-location recording

If you selected multiple locations, it depends on whether you [Select multiple locations to record separate sessions for each location](#) or [Select multiple locations to record a composite session](#) how annotations are stored.

## Separate sessions


The remarks are added to the session that is open on your screen only.

## Composite session


Since all locations have the same Event Log, the remarks are added to all locations.

## procedure

1. Click this button, or press **Ctrl + R**.

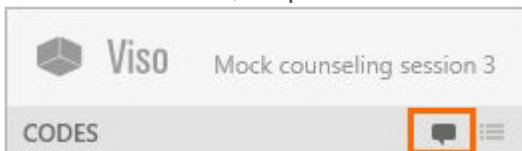

2. Enter the text and click **Add**.

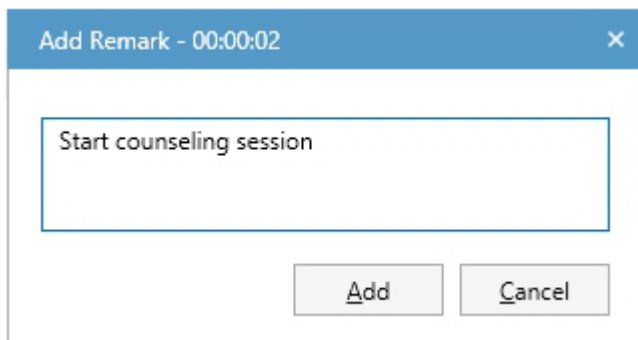

**NOTE** A remark with the location name is added when you switch locations in a composite session.

---

Revision #3

Created 15 January 2024 11:38:49

Updated 11 March 2024 14:26:38