


How Viso annotations appear in The Observer XT

In the Coding Scheme

In The Observer XT, choose **Setup > Open Coding Scheme**, to open the Coding Scheme. Viso subjects appear in the **Subjects** panel. Viso markers are added as Point events in a Start-Stop group. Remarks are not added to the Coding Scheme.

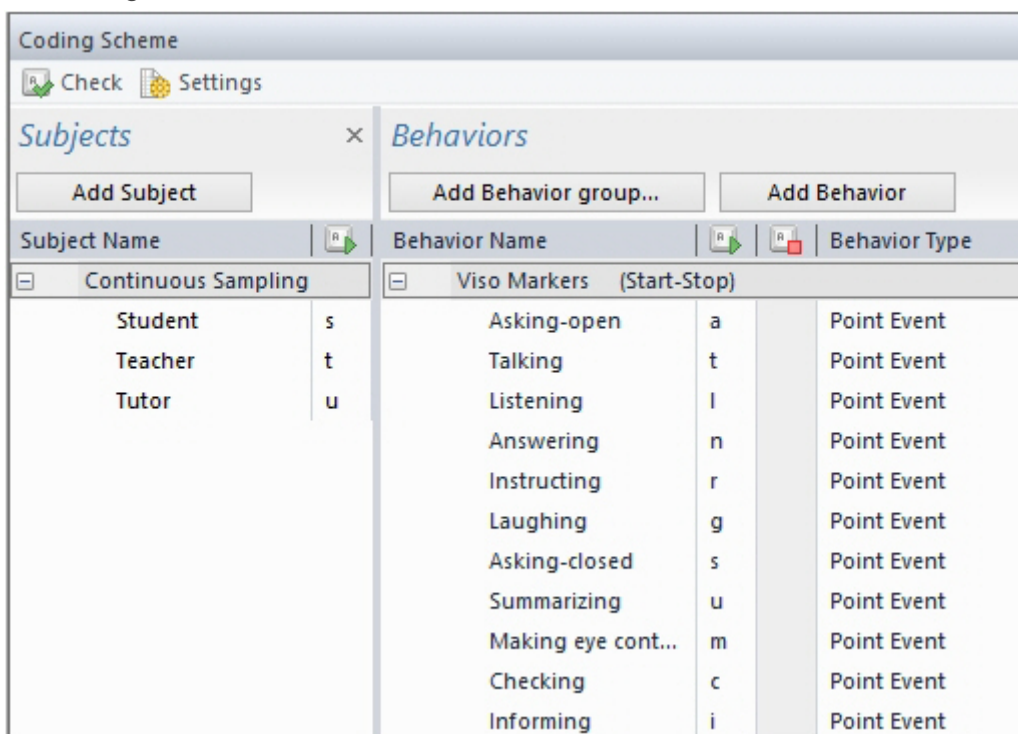

In the observation

Choose **Observe > Observation > Open > [session name and time]** to view the annotations in the Event Log. Viso Subjects are also Subjects in The Observer XT. Markers are point events and are indicated with blue dots. Remarks are imported as comments in the Event Log.

Time	Subject	Behavior	Comment
0.00	Start		
12.06	Student	● Talking	
17.46	Teacher	● Asking-open	
24.34	Student	● Listening	
30.62	Student	● Answering	
39.22	Teacher	● Instructing	
49.82	Student	● Laughing	
55.75	Teacher	● Laughing	
59.31	Student	● Asking-closed	
67.75	Teacher	● Summarizing	
75.03	Student	● Making eye cont	
80.47	Teacher	● Making eye cont	
86.27			Disturbance in back of classroom
99.03	Teacher	● Checking	
106.08	Student	● Summarizing	
108.24	Teacher	● Asking-open	
116.44	Student	● Answering	
125.74	Stop		

To jump to the video frame at which a certain marker or remark was scored, click its row in the Event Log.

Revision #1

Created 15 January 2024 11:38:52

Updated 15 January 2024 11:43:02