

# Introduction

Viso® is a powerful tool to record video and audio in multiple independent locations. Record with up to four Pan-Tilt-Zoom cameras in each room. Control session recording from any computer in your local network, so that the persons under study are not disturbed by the recordings. Annotate the behaviors of the observed persons. Play the recorded sessions back for debriefing. And invite others to view the sessions with you.



## Viso modules

Viso is modular, which means that you can make the setup entirely to your needs.


### Base module

The Viso Base Module allows you to record sessions in a number of locations with up to four cameras in each location. The number of locations is determined by your license. User management ensures that session viewing, recording, or management can only be done by those whose user role allows it.

See [User management](#) for an overview of the rights of each user role.

### Annotation Module

With the Annotation Module you can mark events that are of interest in your session and enter free text as remarks. When debriefing the session, you can easily jump to an episode using the scored markers and remarks.



Team skills training 1

⌕ BACK


EVENTS

TIME	SUBJECT	MARKER	
00:00:12	Student	Talking	●
00:00:17	Teacher	Asking-...	■
00:00:24	Student	Listening	◀
00:00:30	Student	Answeri...	▼
00:00:39	Teacher	Instructi...	+
00:00:49	Student	Laughing	
00:00:55	Teacher	Laughing	
00:00:59	Student	Asking-...	▲
00:01:07	Teacher	Summar...	=
00:01:15	Student	Making...	○
00:01:20	Teacher	Making...	○
00:01:26		Disturba...	■
00:01:39	Teacher	Checking	✕
00:01:46	Student	Summar...	=
00:01:48	Teacher	Asking-...	■
00:01:56	Student	Answeri...	▼

CODES

EVENTS

PARTICIPANTS



TIMELINE


00:00 - 00:10

00:00:00 00:01:00

Student

Teacher

No subject

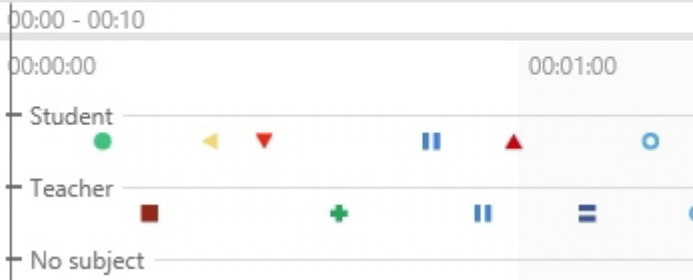

The sessions can be exported to the annotation software The Observer XT, for further analysis.

## Scheduler Module

With this module you can schedule sessions in advance. The Scheduler has several options to customize the view entirely to your needs.

Revision #1

Created 15 January 2024 11:38:43

Updated 15 January 2024 11:43:01