

# Reclaim the role of annotator

## aim

To claim back the role of annotator that you gave to another user, for instance, a colleague or one of your students.

## prerequisites

- Your license includes the Annotation Module.
- You created the session or you received session ownership or you are a Technician or Administrator.

## procedure

1. Open the **Sessions** tab to view the list of all the sessions.
2. Right-click the appropriate session and select **Change annotator**.

NAME	
Interviewing techniques 1	Rename F2
Interviewing techniques 2	Delete Del
Interviewing techniques 3	Change annotator
Test session	Copy videos
Group 1	Import videos
Group 2	Export to The Observer XT
Group 5	Export to CSV

3. Select your name and click **OK**.