

# Review sessions

## aim


To play back recorded sessions. If your license includes the Annotation Module you can also play back the scored markers and events.

## prerequisites

You either have rights to review the session (Technician, Administrator), you created the session (Trainer), you have received ownership or annotator rights or you are invited to view the session.

## procedure


1. Open the **Sessions** tab for a list of all the sessions.

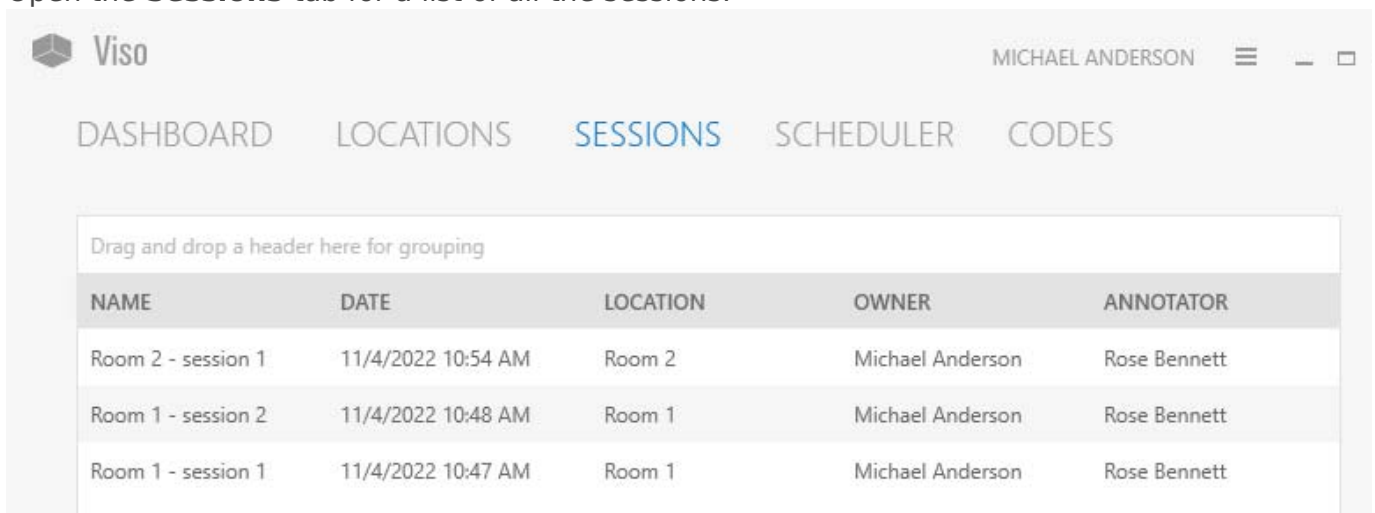

The screenshot shows the Viso application interface. At the top, there is a header bar with the Viso logo on the left and the user name 'MICHAEL ANDERSON' on the right. Below the header, there is a navigation bar with five tabs: 'DASHBOARD', 'LOCATIONS', 'SESSIONS' (which is highlighted in blue), 'SCHEDULER', and 'CODES'. Below the navigation bar, there is a table with the following columns: 'NAME', 'DATE', 'LOCATION', 'OWNER', and 'ANNOTATOR'. The table contains three rows of session data.

NAME	DATE	LOCATION	OWNER	ANNOTATOR
Room 2 - session 1	11/4/2022 10:54 AM	Room 2	Michael Anderson	Rose Bennett
Room 1 - session 2	11/4/2022 10:48 AM	Room 1	Michael Anderson	Rose Bennett
Room 1 - session 1	11/4/2022 10:47 AM	Room 1	Michael Anderson	Rose Bennett

2. Double-click a row to open a session.
3. Click the play button to play it back.

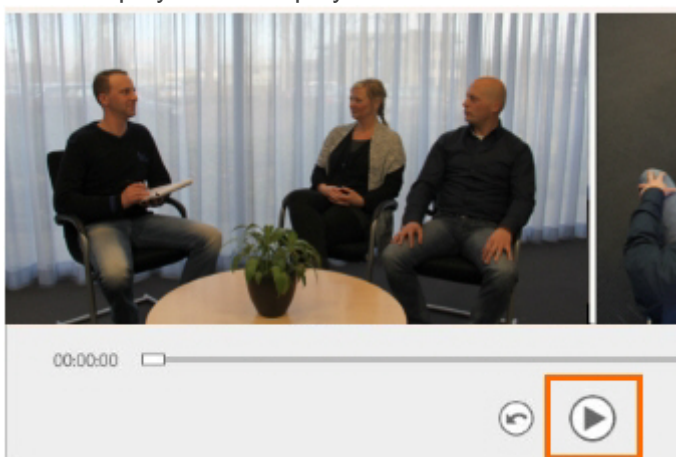

## To review an event

If you want to play back a specific event, set a pre-roll interval to start playing a few seconds prior to the event.

1. Click the menu button, choose **Settings** and then **Playback settings**.
2. Set the duration of the pre-roll interval and click **OK**.

By default the pre-roll interval is set to 5 seconds. The maximum duration you can set is 15 seconds. If you do not want to have a pre-roll, set the value to 0.


## To review a composite session

If your session consists of multiple locations, use the carousel control to view the different locations.

See [To open locations with the carousel control](#)



The Timeline shows when a location change takes place.

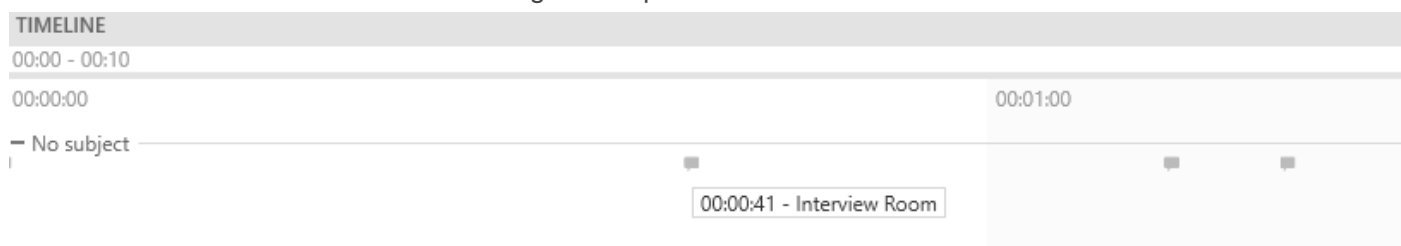

Use the carousel to view the video in the next location, for instance when following a participant from one room to another.

## To change the video layout

Optionally, change the video layout in the same way as with recording sessions.

See [Video layout](#)


## To change the audio

Optionally, change the playback audio in the same way as with recording sessions.


See [Change Audio source and volume in preview](#)

## To review the last time fragment

Click the quick review button to let the video jump back and replay the episode. Click it multiple times to replay a longer episode.

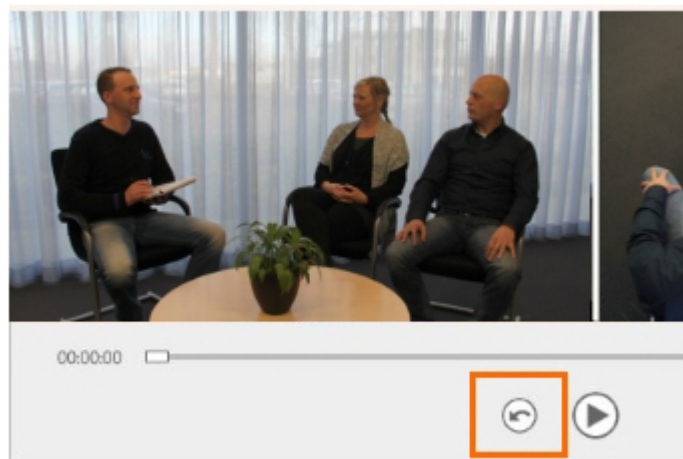

By default, the video jumps back five seconds. To select another time interval, click the **Menu** button at the top of the Viso window and choose **Settings > Playback settings**. Next, specify the interval. You can also control here if you want to pause after quick review or continue the playback.

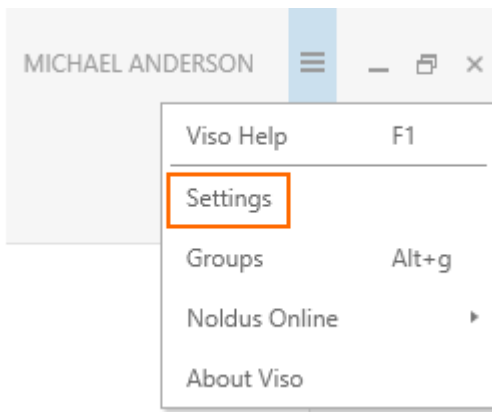

## Session sorting options

Click one of the column headers to sort the sessions, by name, date, location, the person that owns the session or the person that has been assigned as the annotator.

To group sessions, for example by the user that created them, drag that column header (Created by in this example) to the top of the table.

A screenshot of the Viso application interface. The top bar shows the Viso logo and the user's name 'MICHAEL ANDERSON'. Below the top bar is a navigation menu with 'DASHBOARD', 'LOCATIONS', 'SESSIONS' (highlighted in blue), 'SCHEDULER', and 'CODES'. Below the navigation menu is a 'Grouped by:' dropdown menu with 'OWNER' selected and a small upward arrow. Below this is a table with columns: NAME, DATE, LOCATION, OWNER, and ANNOTATOR (with a small upward arrow). The table is grouped by owner, with expandable sections for 'Julius Leurs', 'Michael Anderson', and 'Shahzod Tuychiev'. Each section contains one or more session rows.

NAME	DATE	LOCATION	OWNER	▲ ANNOTATOR
▼ Julius Leurs				
Room 1 - recording 1	9/29/2022 4:22 PM	Room 1	Julius Leurs	Julius Leurs
▼ Michael Anderson				
Room 1 - session 1	11/4/2022 10:47 AM	Room 1	Michael Anderson	Rose Bennett
Room 1 - session 2	11/4/2022 10:48 AM	Room 1	Michael Anderson	Rose Bennett
Room 2 - session 1	11/4/2022 10:54 AM	Room 2	Michael Anderson	Rose Bennett
▼ Shahzod Tuychiev				
Session 1 - room 4	11/3/2022 3:52 PM	Room 4	Shahzod Tuychiev	Shahzod Tuychiev

Unfold a group to see the individual sessions. Click the **red cross** to remove grouping.

A screenshot of the Viso application interface, similar to the previous one, but with the 'Grouped by:' dropdown menu expanded. The dropdown menu shows 'OWNER' with a small red cross icon next to it. Below the dropdown menu is a table with columns: NAME, DATE, and LOCATION.

NAME	DATE	LOCATION
------	------	----------

You can drag multiple column headers to the top of the table to make groups and subgroups of sessions.

---

Revision #6  
Created 15 January 2024 11:38:47  
Updated 31 May 2024 14:16:44 by Niek Wilmink