

Select another Codes list

aim

To choose another Codes list than the one that was used for scoring and annotate with that list.


IMPORTANT To choose another Codes list, you must delete all annotations from the original list first. The deleted annotations cannot be restored.

prerequisites

- Your license includes the Annotation Module.
- You created the session, received session ownership or annotator rights or are a Technician or Administrator.

procedure

1. Open the session.
2. Right-click the markers that were already scored in the Event Log and click **Delete**.
3. Click this button and select another Codes list.

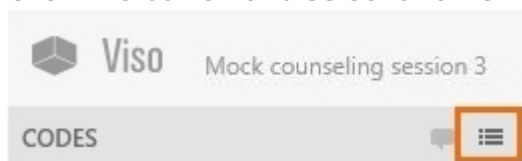

Revision #3

Created 15 January 2024 11:38:50

Updated 8 March 2024 16:36:36