

# Select another Codes list

## aim

To choose another Codes list than the one that was used for scoring and annotate with that list.


**IMPORTANT** To choose another Codes list, you must delete all annotations from the original list first. The deleted annotations cannot be restored.

## prerequisites

- Your license includes the Annotation Module.
- You created the session, received session ownership or annotator rights or are a Technician or Administrator.

## procedure

1. Open the session.
2. Right-click the markers that were already scored in the Event Log and click **Delete**.
3. Click this button and select another Codes list.

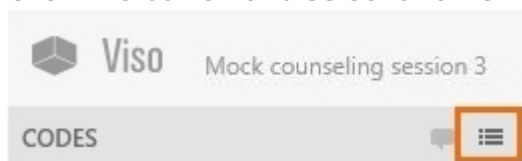

---

Revision #3

Created 15 January 2024 11:38:50

Updated 8 March 2024 16:36:36