Select one location and record a session

aim

To select one location for a single session recording.

prerequisite

Your user role allows session recording (Technician, Administrator, or Trainer).

procedure

1. In the Locations tab, double-click the name of the location in which you want to start recording.

DASHBOARD	LOCATIONS	SESSIONS	CODES		S	
NAME		STATUS				
Control room		Offline				
Demonstration room		Offline	Offline			
Group study room 1.02		Locked	Locked (Rose Bennett)			
Groupb study room 0.101		Online				
Individual study room 1.76		Offline				
Instruction room 0.80		Online				
Instruction room 0.82		Recordi	Recording (Rose Bennett)			
Observation room 1.20		Offline	Offline			
Waiting room		Offline	Offline			

2. A preview of the cameras in the location appears. The toolbar displays the location name.



3. Continue with the third chapter of the Reference (Prepare session recording)

NOTE The location is locked for recording by others at the moment you select it in the Locations list. **NOTE** You can also record a single session with the Viso Web App, see The Viso Web App

Revision #4 Created 15 January 2024 11:38:47 Updated 12 March 2024 09:16:03